



User Manual

OriCell™ Adipose-Derived Stem Cell Growth Medium

Cat. No. GUXMD-90011

PRODUCT DESCRIPTION:

OriCell™ Adipose-Derived Stem Cell Growth Medium consists of optimized Adipose-Derived Stem Cell (ADSC) Basal Medium and pre-selected fetal bovine serum. This product has been developed to optimize the proliferation of adipose-derived stem cells, which can then be expanded through multiple passages while retaining the ability to differentiate into osteoblasts, adipocytes, and chondrocytes.

The product is intended for laboratory research use only. It is not intended for diagnostic, therapeutic, clinical, household, or any other applications.

KIT COMPONENTS:

Adipose-Derived Stem Cell (ADSC) Basal Medium (Cat. No. GUXMD-03011)	440 mL
Adipose-Derived Stem Cell (ADSC)-Qualified Fetal Bovine Serum (Cat. No. GUXMD-05001)	50 mL
Penicillin-Streptomycin	5 mL
Glutamine	5 mL

INSTRUCTIONS:

Preparation of the Complete Medium

1. Prior to use, thaw the ADSC-Qualified Fetal Bovine Serum at 2-8°C overnight or until complete thawed. Gently swirl the bottle to ensure homogeneity. The serum has been heat-inactivated and is ready to use after thawing.



Note: The thawed serum may contain some flocculent precipitates. The presence of these substances in serum does not alter the performance characteristics of the product. It is not recommended to filter the serum alone to remove these precipitates. Doing so may result in the loss of some serum nutrients.

2. About 30 minutes prior to use, thaw the Penicillin-Streptomycin solution and Glutamine solution at room temperature. Gently swirl the vials to ensure homogeneity.
3. Disinfect the external surfaces of the bottles/vials for every component in the kit with 70% v/v ethanol. Allow ethanol to evaporate.
4. Aseptically open the bottles/vials inside a laminar flow hood.
5. Transfer the entire amount of ADSC-Qualified Fetal Bovine Serum, Penicillin-Streptomycin solution, and Glutamine solution into the ADSC Basal Medium.
6. Rinse each vial with a small amount of basal medium. Subsequently transfer the rinse medium back into the bottle of basal medium.
7. Gently swirl the fully supplemented (complete) medium to ensure a homogeneous

mixture. The complete medium is now ready to use.



Note: Although each component in this kit is supplied sterile, it is strongly recommended to filter the fully supplemented (complete) medium.

STABILITY AND STORAGE:

All products should be stored in the dark. Adipose-Derived Stem Cell Basal Medium is stable at 2-8°C for up to one year. Other components are stable at -20°C for up to two years.

These products should be discarded beyond the labeled expiration date. Once prepared, the fully supplemented (complete) medium can be stored for up to one month when stored in the dark at 2-8°C.

For optimal performance, repeated warm-cooling and freeze-thawing should be avoided.

QUALITY CONTROL:

OriCell™ Adipose-Derived Stem Cell Growth Medium has been tested for performance on adipose-derived stem cells. The standard evaluation includes:

- Sterility test (bacteria, fungi, and mycoplasma)
- pH test
- Osmolality
- Endotoxin

RELATED PRODUCTS:

Product	Catalog Number
OriCell™ Wistar Rat Adipose-Derived Mesenchymal Stem Cells	RAWMD-01001
OriCell™ SD Rat Adipose-Derived Mesenchymal Stem Cells	RASMD-01001
OriCell™ Rabbit Adipose-Derived Mesenchymal Stem Cells	RBXMD-01001
OriCell™ Dog Adipose-Derived Mesenchymal Stem Cells	CAXMD-01001

Cyagen Biosciences reserves all rights on the technical documents of its OriCell™ cell culture products. No part of this document may be reproduced or adapted for other

purposes without written permission from Cyagen Biosciences.